Caramuel 38, 28011 Madrid • Tel. 91 464 94 50 - 91 464 36 00
Fax. 91 464 62 58 • www.f-soria.es

FICHA TÉCNICA: 770646

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REFERENCIAS BIBLIOGRAFICAS

1. Baird-Parker, A.C. 1962. An improved diagnostic and selective medium for isolating coagulase-positive staphylococci.. J. Appl. Bacteriol. 25: 12-19.
2. Lancette, G.A., and R.W. Bennett. 2001. *Staphylococcus aureus* and staphylococcal enterotoxins. In: Downes, F.P., and K. Ito (ed.). Compendium of methods for the microbiological examination of foods, 4th edition. American Public Health Association (APHA). Washington, D.C. USA.
3. Flowers, R.S., W. Andrews, C.W. Donnelly, and E. Koenig. 1993. Pathogens in milk and milk products, p. 103-212. In: R.T. Marshall (ed.). Standard methods for the examination of dairy products, 16th ed. American Public Health Association, Washington DC, USA.

4. Association of Official Analytical Chemists. 1995. Bacteriological Analytical Manual. 8th edition. AOAC International. Arlington, VA, USA.
5. United States Pharmacopeial Convention, Inc. 1995. The United States Pharmacopeia, 23rd edition. The United States Pharmacopeial Convention, Inc. Rockville, MD, USA.
6. Eaton, A.D., L.S. Clesceri, and A.E. Greenberg (ed.). 1995. Recreational waters, p. 9.26 – 9.27. *In:* Standard methods for the examination of water and wastewater, 19th edition. American Public Health Association, Washington DC, USA.
7. MacFaddin, J.F. 1985. Media for the isolation – cultivation – maintenance of medical bacteria. Volume 1. Williams and Wilkins, Baltimore, London
8. Chapin, K.C., and T.-L. Lauderdale. 2003. Reagents, stains, and media: bacteriology. *In:* Murray, P. R., E. J. Baron, J.H. Jorgensen, M. A. Pfaller, and R. H. Tenover (ed.). Manual of clinical microbiology, 8th ed. American Society for Microbiology, Washington, D.C.
9. Bannerman, T.L. 2003. *Staphylococcus*, *Micrococcus*, and other catalase-positive cocci that grow aerobically. *In:* Murray, P. R., E. J. Baron, J.H. Jorgensen, M. A. Pfaller, and R. H. Tenover (ed.). Manual of clinical microbiology, 8th ed. American Society for Microbiology, Washington, D.C.

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